

Inert Draw-out plate Lens

For use with 5973 Inert MS sources.

5973 Inert Draw-out plate Lens: For use with 5973 Inert MS Source only.

Note: How can you tell if you have an inert source? There are several ways to identify an inert source. First, the screws that secure the inert source (repeller, filaments and lens set screw) are gold. There is a spacer/washer between the repeller and source body. Finally, the filaments are different and the affected parts (source body, repeller assembly and draw out plate) are marked by a dimple.

Warning: It is very important NOT to mix and match parts from a standard source, extended linearity source and the inert source or problems with unpredictable results can occur.

Part Number	Part ID
G2589-20100	Draw-out plate lens

Associated Inert MS source parts:

G2590-60053	High temperature Filaments
G2589-60102	Complete Repeller Assembly (Includes heater & Sensor)
G1999-20021	Gold Source screws (For securing Repeller and Filaments to Source)
G1999-20022	Gold lens stack set screw

This document is believed to be accurate and up-to-date. However, Agilent Technologies, Inc. cannot assume responsibility for the use of this material. The information contained herein is intended for use by informed individuals who can and must determine its fitness for their purpose.