




Agilent Automation Application Highlight

Promega Wizard SV96 PCR cleanup on the BenchCel Workstation

PCR Clean-up Workstation

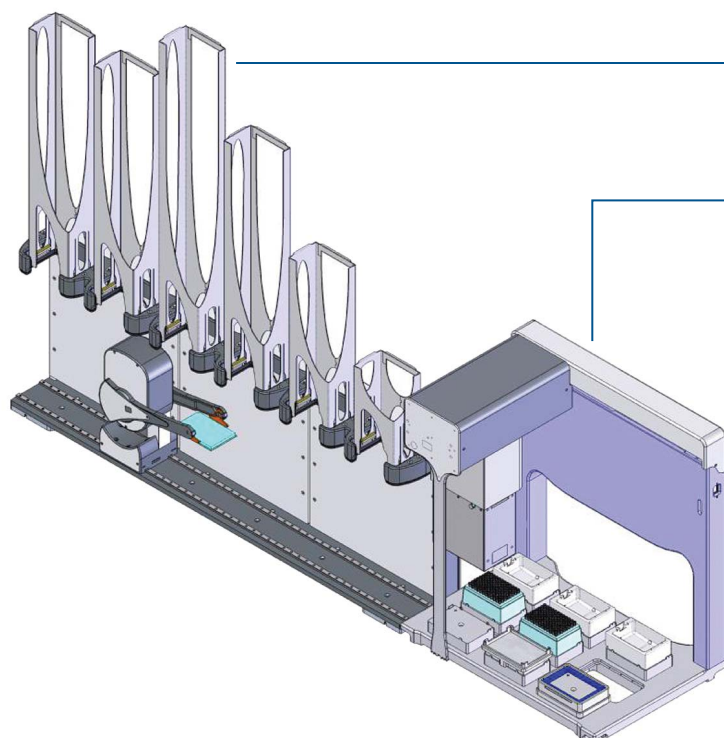

Agilent BenchCel Microplate Handler

Storage of sample plates, Promega Wizard SV96 Binding and Elution plates, and tipboxes.

Agilent Bravo Liquid Handler

Addition of sample, binding buffer, wash solutions, and elution buffer to binding and elution plates. On-deck Bravo Vacuum Filtration Station for sample washes and elution.

A Benchtop Solution:

Agilent Automation provides a complete benchtop solution for PCR Clean-up.

The combination of the Bravo Automated Liquid Handling Platform and the BenchCel Microplate Handler allows for a flexible, efficient and compact approach to the automation of PCR Clean-up. The BenchCel Microplate Handler has capacity to deliver 15 kits to the Bravo Liquid Handler for processing. The Bravo Liquid Handler automatically assembles and disassembles the vacuum manifold through multiple elution cycles without user interaction for true walk-away performance.


Agilent Automation Benefits:

- **Plate Handling:** Use most 1536-, 384-, and 96-well microplates, PCR plates, tube racks, and lids.
- **Throughput:** A PCR plate is processed every 20 minutes.
- **Scalability:** Additional stacks or BenchCels can increase capacity and throughput as needed.
- **Capability:** The Bravo automatically assembles and disassembles the vacuum manifold.
- **Intelligent:** VWorks Automation Control software schedules protocol tasks, incubation steps, and reagent order-of-addition for maximum throughput and walk-away time.



Agilent Technologies

Overview of the Promega Wizard SV96 PCR Clean-up Assay

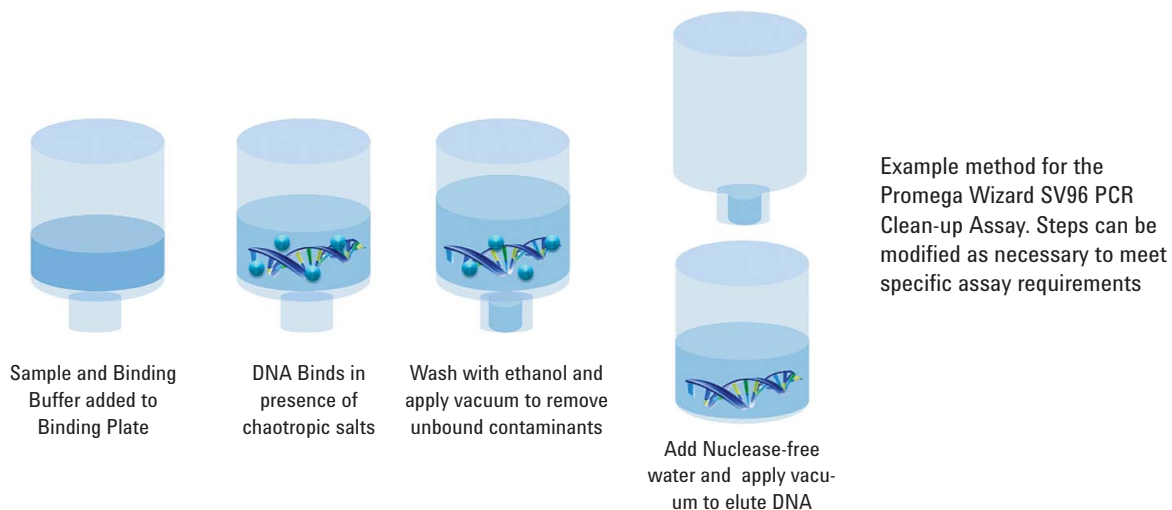

Configuration Options

Agilent BenchCel Workstations offer the flexibility of adding and reconfiguring options to meet your workflow requirements.

- **Integration:** Add a PlateLoc Microplate Sealer, Microplate Labeler, Microplate Centrifuge, or a third-party device to the free end of the BenchCel Microplate Handler.
- **Accessories:** The Bravo Automated Liquid Handling Platform has a growing list of interchangeable deck accessories. Add a Barcode Reading Station for plate tracking or an Orbital Shaking station for efficient mixing.
- **More:** Speak with one of our product specialists or application specialists to determine which configuration will work best for you.

Additional Services

All BenchCel Workstations include installation and product familiarization along with options to purchase additional services to meet your specific needs.

These options include:

- **Consulting:** Work with a team of experts to determine the best solution for your laboratory.
- **Application Development:** Let our application specialists develop and test system protocols for your applications.
- **More:** Speak with one of our product specialists regarding any additional support you require to ensure our solution meets all of your needs.

www.agilent.com/lifesciences/automation

Agilent Technologies shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2009
Published in the USA
October 26, 2009
5990-4951EN



Agilent Technologies