

G2350A Emission source assembly

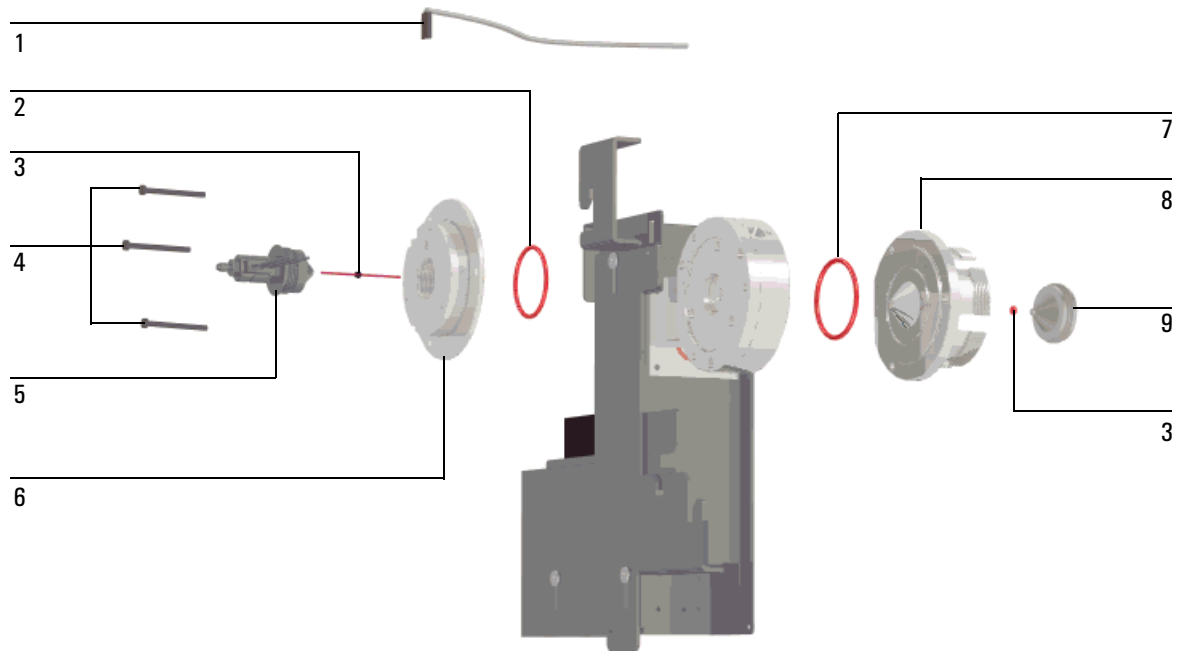


Figure 157. Emission source assembly, part no. G2350-60560

Number	Description	Part number	Qty/Assy
	Emission source assembly	G2350-60560	–
	RP-emission source assembly	G2350-60900	–
1	Cavity heater/PRT assembly	G2350-60930	1
2	GC chamber O-ring	0905-1159	1
3	Discharge tube kit	G2350-61090	1
4	RP-MS M5×3.80 35mm	0515-0110	3
5	Gas union assembly	G2350-61190	1
6	RP-heater chamber assembly	G2350-60190	1
7	Optical side water chamber O-ring	0905-1160	1
8	Optical side water chamber	G2350-20520	1
9	Exit chamber	05921-22240	1
NS	Belleville washers	3050-1907	4
NS	Set screw	0515-2432	